
bmlab Documentation

Release v0.8.0.post8

Paul Müller

Apr 17, 2023

CONTENTS:

1	Getting started	3
1.1	Installation	3
1.2	Basic usage	3
1.3	Citing bmlab	4
2	Code reference	5
3	Development	7
3.1	Tests	7
3.2	Making a new release	7
3.3	Continuous integration	7
4	Changelog	9
4.1	version ## 0.7.0 - 2022-09-28	11
4.2	version	11
4.3	version ### Added	11
4.4	version - Implement 4-peak fitting #167	11
4.5	version	11
4.6	version ## 0.6.1 - 2022-09-16	11
4.7	version	11
4.8	version ### Fixed	11
4.9	version - Fix get_frequencies_by_time #165	11
4.10	version - Fix evaluating files without calibration #166	11
4.11	version	11
4.12	version ## 0.6.0 - 2022-09-14	11
4.13	version	11
4.14	version ### Added	11
4.15	version - Implement exporting 3D data #162	11
4.16	version	11
4.17	version ### Changed	11
4.18	version - Fit Brillouin spectrum on frequency axis #158	11
4.19	version	11
4.20	version ### Fixed	11
4.21	version - Fix parameter label in csv files #163	11
4.22	version	11
4.23	version ## 0.5.1 - 2022-09-02	11
4.24	version	11
4.25	version ### Fixed	11
4.26	version - Correctly initialize setup on new file load #156	11
4.27	version	11

4.28	version ## 0.5.0 - 2022-09-02	11
4.29	version	11
4.30	version ### Added	11
4.31	version - Also store Rayleigh peak shift #153	11
4.32	version - Write scale bar to Brillouin Tiff file #154 #155	11
4.33	version	11
4.34	version ## 0.4.0 - 2022-08-29	11
4.35	version	11
4.36	version ### Added	11
4.37	version - Allow to specify calibration temperature #143	11
4.38	version - Export all peaks from multi-peak fits #146	11
4.39	version - Also store peak fit offsets #150 #151	11
4.40	version - Allow to get fits #152	11
4.41	version	11
4.42	version ### Changed	11
4.43	version - Update scipy #142	11
4.44	version	11
4.45	version ### Fixed	11
4.46	version - Handle indices as int when getting key #144	11
4.47	version - Fix exporting fluorescence w/o scale calibration #145	11
4.48	version - Don't overwrite bounds when setting peak number #148	11
4.49	version - Don't fail to set regions w/o calibration #149	11
4.50	version	11
4.51	version ## 0.3.0 - 2022-08-17	11
4.52	version	11
4.53	version ### Added	11
4.54	version - Allow to export all Brillouin parameters acquired #141	11
4.55	version	11
4.56	version ### Fixed	11
4.57	version - Align export parameter names with rest of library #140	11
4.58	version	11
4.59	version ## 0.2.4 - 2022-08-12	11
4.60	version	11
4.61	version ### Fixed	11
4.62	version - Don't fail on empty arc #138	11
4.63	version - Fix issues with flake8 v5, lock scipy version #139	11
4.64	version	11
4.65	version ## 0.2.3 - 2022-07-01	11
4.66	version	11
4.67	version ### Fixed	11
4.68	version - Fix exporting on macOS #135	11
4.69	version - Resolve paths to ensure proper folder handling #137	11
4.70	version - Minor cleanup of test_serializer.py #136	11
4.71	version	11
4.72	version ## 0.2.2 - 2022-06-29	11
4.73	version	11
4.74	version ### Fixed	11
4.75	version - Explicitly import pdfpages to solve pyinstaller issue #134	11
4.76	version	11
4.77	version ## 0.2.1 - 2022-06-28	11
4.78	version	11
4.79	version Re-release to trigger PyPi upload. No changes to 0.2.0.	11
4.80	version	11
4.81	version ## 0.2.0 - 2022-06-27	11

4.82	version	11
4.83	version ### Added	11
4.84	version - Implement exporting Brillouin and fluorescence maps as images #116	11
4.85	version - Implement exporting Brillouin data as CSV file #131	11
4.86	version - Implement retrieving fluorescence images #115	11
4.87	version - Compensate drift by shifting evaluate regions #114	11
4.88	version - Allow to get image binning #123 and binning factor #124	11
4.89	version	11
4.90	version ### Changed	11
4.91	version - Evaluate in the temporal order of acquisition #113	11
4.92	version - Reduce code duplication #112	11
4.93	version - Optimize calculation of get_arc_from_circle_phis #125	11
4.94	version - Account for binning when extracting #127	11
4.95	version - Allow to set all points at once when extracting #129	11
4.96	version - Only update the positions for the given key #130	11
4.97	version	11
4.98	version ### Fixed	11
4.99	version - Improve the spectra extraction #120 #132	11
4.100	version - Ensure discretize_arc returns arc within image #121	11
4.101	version - Only use linear frequency interpolation #122	11
4.102	version - Don't crash finding all points on empty calibrations #128	11
4.103	version - Don't crash evaluation when no spectra is returned #133	11
4.104	version	11
4.105	version ## 0.1.10 - 2022-05-12	11
4.106	version	11
4.107	version ### Changed	11
4.108	version - Raise proper exceptions if source or session files are lacking #106 #111	11
4.109	version - Improve exact circle calculation #108	11
4.110	version	11
4.111	version ### Fixed	11
4.112	version - Don't fail hard for invalid circle fits #107	11
4.113	version - Don't crash when changing number peaks during evaluation #109	11
4.114	version - Don't fail hard for aborted measurements #110	11
4.115	version	11
4.116	version ## 0.1.9 - 2022-05-06	11
4.117	version	11
4.118	version ### Fixed	11
4.119	version - Don't fail hard for values outside interpolation range #104	11
4.120	version - Fix calculating and handling bounds #105	11
4.121	version	11
4.122	version ## 0.1.8 - 2022-05-03	11
4.123	version	11
4.124	version ### Added	11
4.125	version - Support multi-peak fitting for measurement data #99 #102	11
4.126	version - Adjust evaluation model to store multi-peak fit results #96	11
4.127	version - Implement getting correct multi-peak fit data #97	11
4.128	version - Add an example script for exporting 2D data #101	11
4.129	version	11
4.130	version ### Changed	11
4.131	version - Interpolate to find index for given frequency #100	11
4.132	version - Improve the calibration peak finding #103	11
4.133	version	11
4.134	version ## 0.1.7 - 2022-04-22	11
4.135	version	11

4.136 version ### Added	11
4.137 version - Allow to set calibration frequencies #87	11
4.138 version - Add functions to get image count #92	11
4.139 version - Allow to sort calibration keys by time #93	11
4.140 version - Add function to get expected calibration frequencies #95	11
4.141 version	11
4.142 version ### Changed	11
4.143 version - Adjust required peak height to signal level #86	11
4.144 version	11
4.145 version ### Fixed	11
4.146 version - Delete VIPA params when clearing calibration #88	11
4.147 version - Fix issue when only one calibration sample is used #89	11
4.148 version - Only clear correct fit, add test for FitSet #94	11
4.149 version	11
4.150 version ## 0.1.6 - 2022-04-11	11
4.151 version	11
4.152 version ### Fixed	11
4.153 version - Better handle broken eval data files #85	11
4.154 version	11
4.155 version ## 0.1.5 - 2022-04-11	11
4.156 version	11
4.157 version ### Fixed	11
4.158 version - Don't fail on empty repetitions #84	11
4.159 version	11
4.160 version ## 0.1.4 - 2022-03-22	11
4.161 version	11
4.162 version ### Fixed	11
4.163 version - Bring back multiprocessing support #82	11
4.164 version	11
4.165 version ## 0.1.3 - 2022-03-21	11
4.166 version	11
4.167 version ### Fixed	11
4.168 version - Improve circle fit for points on straight line #80	11
4.169 version - Remove multiprocessing support for now #81	11
4.170 version	11
4.171 version ## 0.1.2 - 2022-03-15	11
4.172 version	11
4.173 version ### Fixed	11
4.174 version - Maintain order of parameter keys #78	11
4.175 version	11
4.176 version ## 0.1.1 - 2022-03-15	11
4.177 version	11
4.178 version ### Added	11
4.179 version - Provide session files for source files #74	11
4.180 version	11
4.181 version ### Fixed	11
4.182 version - Ensure we always get the expected data shape #75	11
4.183 version	11
4.184 version ## 0.1.0 - 2022-03-14	11
4.185 version	11
4.186 version ### Added	11
4.187 version - Implement getting results maps #70	11
4.188 version - Save bmlab version in session file #71	11
4.189 version - Allow to open session files directly #72	11

4.190 version	11
4.191 version ## 0.0.13 - 2022-02-21	11
4.192 version	11
4.193 version ### Fixed	11
4.194 version - Fix lorentz fit for special case #68	11
4.195 version	11
4.196 version ## 0.0.12 - 2022-02-16	11
4.197 version	11
4.198 version ### Added	11
4.199 version - Add example for standalone evaluation #61	11
4.200 version - Allow finding peaks in extraction view for all calibrations #65	11
4.201 version - Allow to rotate the image by negative number #67	11
4.202 version	11
4.203 version ### Changed	11
4.204 version - Improve the peak finding for calibration spectra #63	11
4.205 version - Code cleanup #66	11
4.206 version	11
4.207 version ### Fixed	11
4.208 version - Correctly close and join mp pool #60	11
4.209 version - Fix switched Rayleigh and Brillouin regions in test #62	11
4.210 version	11
4.211 version ## 0.0.11 - 2022-02-01	11
4.212 version	11
4.213 version ### Added	11
4.214 version - Implement peak finding for calibration spectra #56	11
4.215 version - Allow to add regions by frequency #59	11
4.216 version	11
4.217 version ### Changed	11
4.218 version - Don't show runtime warning for fitting #54	11
4.219 version - Calculate derived values during evaluation #55	11
4.220 version - Bundle auto evaluation in a general controller #58	11
4.221 version - Code cleanup #52 #57	11
4.222 version	11
4.223 version ### Fixed	11
4.224 version - Correct bmlab citation hint #51	11
4.225 version - Don't fail when not enough peaks for frequency axis #53	11
4.226 version	11
4.227 version ## 0.0.4	11
4.228 version	11
4.229 version - setup: use “pytest” command instead of deprecated “setup.py test”	11
4.230 version - get metadata from HDF file	11
4.231 version - set image orientation	11
4.232 version	11
4.233 version ## 0.0.3	11
4.234 version	11
4.235 version - testing connection with BMicro	11
4.236 version	11
4.237 version ## 0.0.2	11
4.238 version	11
4.239 version - CI test release	11
4.240 version	11
4.241 version ## 0.0.1	11
4.242 version	11
4.243 version - dummy release	11

4.244 version	11
5 Bibliography	13
6 Indices and tables	15
Python Module Index	17
Index	19

This is bmlab, a Python library for Brillouin microscopy data analysis. This is the documentation of bmlab version v0.8.0.post8.

GETTING STARTED

1.1 Installation

To install bmlab, use one of the following methods:

- from PyPI: pip install bmlab
- from sources: pip install .

1.2 Basic usage

```
import pathlib

from bmlab.controllers import Controller
from bmlab.models import Orientation
from bmlab.models.setup import AVAILABLE_SETUPS

filepath = pathlib.Path(__file__).parent.parent / 'tests' / 'data' / 'Water.h5'
setup = AVAILABLE_SETUPS[0]
orientation = Orientation(rotation=1, reflection={
    'vertically': False, 'horizontally': False
})

# Frequency ranges to evaluate in Hz
brillouin_regions = [(4.0e9, 6.0e9), (9.0e9, 11.0e9)]
rayleigh_regions = [(-2.0e9, 2.0e9), (13.0e9, 17.0e9)]

# The number of Brillouin peaks to fit. Can be set to 2
# for samples where contributions to the Brillouin shift
# from different materials are expected in the same focal
# volume, e.g. droplets.
nr_brillouin_peaks = 1

# Bounds for the multi-peak fit, only considered for nr_brillouin_peaks = 2.
# Possible options: 'min', 'max', '-inf', 'inf' or number in GHz.
multi_peak_bounds = [
    ['min', 5.0],
    [5.5, 'inf']
]
```

(continues on next page)

(continued from previous page)

```
# This condition is necessary due to
# bmlab using multiprocessing!
if __name__ == '__main__':
    session = Controller().evaluate(
        filepath,
        setup,
        orientation,
        brillouin_regions,
        rayleigh_regions,
        nr_brillouin_peaks,
        multi_peak_bounds
    )
    session.save()
```

1.3 Citing bmlab

If you use bmlab in a scientific publication, please cite it with:

bmlab developers (2022), bmlab version X.X.X: Python library for the post-measurement analysis of Brillouin microscopy data sets [Software]. Available at <https://github.com/BrillouinMicroscopy/bmlab>.

CODE REFERENCE

```
class bmlab.file.BrillouinFile(path)
```

Load a HDF file with Brillouin microscopy data.

Parameters `path` (*Path*) – path of the file to load

Raises `OSError` – when trying to open non-existing or bad file

```
get_repetition(repetition_key, mode='Brillouin')
```

Get a repetition from the data file based on given key.

Parameters

- `repetition_key` (*str*) – key to identify the repetition in the Brillouin group
- `mode` (*str*) – the mode to look at

Returns `out` – the repetition

Return type *Repetition*

```
repetition_count(mode='Brillouin')
```

Get the number of repetitions in the data file.

Returns `out` – Number of repetitions in the data file

Return type *int*

```
repetition_keys(mode='Brillouin')
```

Returns list of keys for the various repetitions in the file.

Returns `out`

Return type list of str

```
class bmlab.file.Repetition(repetition_group, file)
```

Creates a repetition from the corresponding group of a HDF file.

Parameters `repetition_group` (*HDF group*) – The HDF group representing a Repetition. Consists of payload, calibration and background.

```
class bmlab.file.Payload(payload_group, repetition)
```

Creates a payload representation from the corresponding group of a HDF file.

Parameters `payload_group` (*HDF group*) – The payload of a repetition, basically a set of images

```
class bmlab.file.Calibration(payload_group, repetition)
```

Creates a calibration representation from the corresponding group of a HDF file.

Parameters `payload_group` (*HDF group*) – Calibration data of a repetition from an HDF file.

Module for interacting with files containing Brillouin microscopy data.

NOTE: Ideally, the users of bmlab should not have to know about the file format in which the data are stored. So, if possible, do not expose HDF objects to the outside (like BMicro).

exception `bmlab.file.BadFileException`

DEVELOPMENT

For the general workflow, please refer to the [BMicro docs](#).

3.1 Tests

We try to adhere to test-driven development. Please always write test functions for your code. Please make sure the `pytest` package is installed:

```
pip install pytest
```

You can run all tests via:

```
py.test tests
```

3.2 Making a new release

The release process of bmlab is completely automated. All you need to know is that you have to create an incremental tag on the main branch:

```
git tag -a "0.1.3"  
# or (if you have set up PGP)  
git tag -s "0.1.3"  
# and finally  
git push --tags
```

For more information on how automatic deployment to PyPI works, please read on.

3.3 Continuous integration

The following things are automated:

- `pytest` and `flake8` on Linux, macOS, and Windows via GitHub Actions: <https://github.com/BrillouinMicroscopy/bmlab/actions?query=workflow%3AChecks>

You should always check that all checks pass before you merge a pull request (A green state on your local machine does not mean a global green state).

- automatic deployment to PyPI on tag creation via GitHub Actions: <https://github.com/BrillouinMicroscopy/bmlab/actions?query=workflow%3A%22Release+to+PyPI%22>

Paul Müller created the `bmlab` package on PyPI and gave the user `ci_bm` permission to upload new releases. The password for this user is an organization secret.

- documentation is built automatically (for all tags and for the latest commit to the main branch) on readthedocs: <https://readthedocs.org/projects/bmlab/builds/>
- coverage statistics are done with codecov: <https://codecov.io/gh/BrillouinMicroscopy/bmlab>

Please try stay above 90% coverage.

Badges for all of these CI tasks are in the main `README.rst` file.

**CHAPTER
FOUR**

CHANGELOG

List of changes in-between bmlab releases.

4.1 version ## 0.7.0 - 2022-09-28

4.2 version

4.3 version ### Added

4.4 version - Implement 4-peak fitting #167

4.5 version

4.6 version ## 0.6.1 - 2022-09-16

4.7 version

4.8 version ### Fixed

4.9 version - Fix get_frequencies_by_time #165

4.10 version - Fix evaluating files without calibration #166

4.11 version

4.12 version ## 0.6.0 - 2022-09-14

4.13 version

4.14 version ### Added

4.15 version - Implement exporting 3D data #162

4.16 version

4.17 version ### Changed

4.18 version - Fit Brillouin spectrum on frequency axis #158

4.19 version

4.20 version ### Fixed

4.21 version ## 0.7.0 - 2022-09-28

4.21 version - Fix parameter label in csv files #163

4.22 version

**CHAPTER
FIVE**

BILBLIOGRAPHY

**CHAPTER
SIX**

INDICES AND TABLES

- genindex
- modindex
- search

PYTHON MODULE INDEX

b

bmlab.file, 6

INDEX

B

`BadFileException`, 6
`bmlab.file`
 `module`, 6
`BrillouinFile` (*class in bmlab.file*), 5

C

`Calibration` (*class in bmlab.file*), 5

G

`get_repetition()` (*bmlab.file.BrillouinFile method*), 5

M

`module`
 `bmlab.file`, 6

P

`Payload` (*class in bmlab.file*), 5

R

`Repetition` (*class in bmlab.file*), 5
`repetition_count()` (*bmlab.file.BrillouinFile method*), 5
`repetition_keys()` (*bmlab.file.BrillouinFile method*),
 5